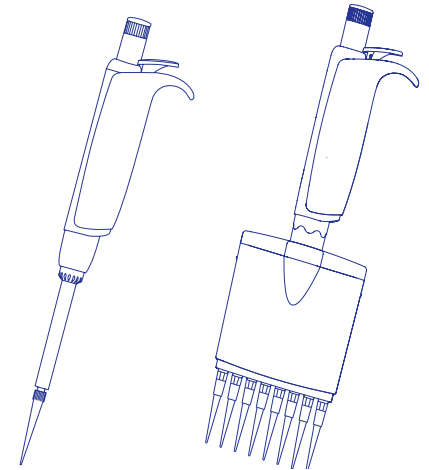
ACURA® manual 815 / 825 / 835 / 855

DIGITAL READING PIPETTES PIPETTEN MIT DIGITAL-DISPLAY PIPETTES À LECTURE DIGITALE

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⚠ SAFETY PRECAUTIONS

- Read operating instructions carefully; keep available for future reference.
- Observe manufacturer's instructions and guidelines for operation and maintenance.
- Refer to and follow regulations about handling of potentially hazardous reagents.
- Before use, check instrument for good working conditions.
- Do not use harsh chemicals (such as acetone) to clean instrument.
- This product should be used only for its intended purpose.
- Mind possible hand fatigue during serial pipetting and its medical consequences (such as repetitive strain injuries RSI).

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DESCRIPTION (Fig. 1)

The Acura® manual is an air displacement micropipette line intended for the accurate, precise and easy pipetting. The single channel models cover volumes ranging from 0.1 µl to 10 mL. Available in ten models with 8 or 12 channels, the multichannels are perfectly adapted for a continuous microplate pipetting in the 0.5 to 350 µL range. Materials have been carefully selected to ensure durability and optimal resistance to aggressive media. The unit is autoclavable at 121°C fully assembled. Major features include digital display window (E) for permanent volume visibility and integrated, adjustable tip ejector Justip™ (H) activated by separate push button (C). Ergonomics and pipette activation provide for an ultra-soft operation. Calibration easily performed by laboratory personal with *swift-set* system (F).

OPERATION (Fig. 1)

Holding the pipette

Place finger rest (D) on the phalanx of forefinger. Ease of access of both plunger button (B) and ejector push button (C) allow natural and comfortable thumb movement. The multichannel casing (K) revolves to allow the selection of the best working position.

Volume setting

Set desired volume of variable models by rotating push button (B) clockwise to decrease the volume and vice versa, until it appears in display window (E). Set volume is exactly indicated by the top of the pointer. Precise click-stops of the micrometric mechanism and a free rotating *smartie* colour cap (A) on top of button prevent any volume alteration during pipetting.

Note: If the letter E appears beside digits in the display window, the set volume is no longer within working range of the pipette. Overturning the micrometric screw may impair the volume setting mechanism.

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Tip selection

Disposable Qualitix® – with or without filter – were carefully designed to provide the best pipette and tip combination for all types of applications. Regular tips are autoclavable at 121°C. Filter tips are either autoclavable or supplied in a sterile version. Use tips supplied by manufacturer, or similar (see tip selection on p II). Detailed flyer available on request.

Apply new tip as often as deemed necessary in regard with change of sample nature, droplet accumulation after blow out, or foam formation. When using tip racks, place the instrument on tips, press firmly to ensure a good tightness. This is particularly important for multichannel models.

Note: Enhance precision by filling new tip at least once before sampling.

Pasteur pipette (835 model only)

Glass Pasteur pipettes are of advantage in handling PP affecting solvents. Optional adapter-nozzles (cat. No. 1.835.631, 2 mL or 1.835.633, 5 mL), accommodate standard 2 mL Pasteur pipettes (ext. Ø 6.5-7.2 mm) in addition to Polypropylene tips. Refer to instructions for disassembling volumetric module for changing the adapter nozzle. Performance shown on QC certificate are obtained with PP tips, recalibrate if necessary with Pasteur pipette.

Note: Do not set volume higher than 2 mL.

Tip ejection (Fig. 2 and 3)

The patented Justip™ ejection system* allows instant height adjustment (± 2 mm) of ejector for best fitting of tip used (optimal distance). Set correct position (~ 0.5 mm space between tip and ejector) by rotating ejector screw ($< LO - HI >$). Click stops prevent any unwanted change while pipetting.

A special design ensure sequential ejection on multichannel models.

Protection filter (835 models)

Macropipettes (2, 5 and 10 mL models) accommodate a nozzle protection filter, which drastically reduces risks of barrel and plunger contamination. Its use is particularly recommended with large volumes and/or when handling biohazardous, radioactive or highly aggressive liquids. Change filter if wet or contaminated. Filter must not be autoclaved.

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*Socorex patented 6, 833, 114B1 (US Patent)

PIPETTING OPERATION

Before pipetting, tightly fit new and clean tips on the nozzle.

Forward pipetting (Fig. 4)

Exact volume desired is aspirated, then dispensed.

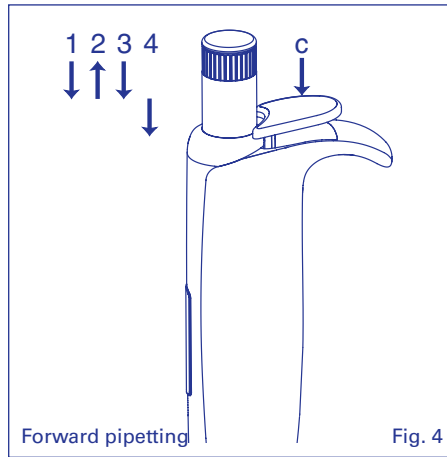
Phase 1 Press plunger button to intermediate stop ending calibrated stroke.

Phase 2 Immerse tips vertically (2 - 3 mm) into liquid, wait 2 sec. and smoothly release plunger button. Withdraw tips from liquid, without touching container wall.

Phase 3 Place tips against wall of receiving vessel, press plunger button to first stop and let liquid out.

Phase 4 Press button to second stop; blow out will expel remaining droplets. Withdraw tips from receiving vessel, sliding it over 10-15 mm along container wall.

Tip ejection Remove used tips by pressing ejector button (C). Change tips as required.

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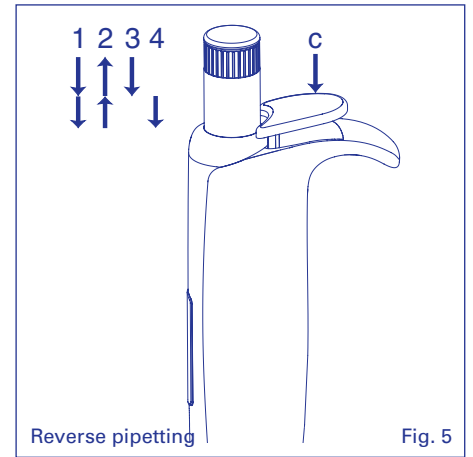
Reverse pipetting (Fig. 5)

Liquid is aspirated in excess of the selected volume but only the set volume is dispensed.

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Phase 3 Place tips against wall of receiving vessel, depress plunger button to first stop only and let liquid out.

Phase 4 Withdraw tips from receiving vessel, sliding it over 10 -15 mm along container wall and start operation to phase 2 again, or discard residual liquid by Pressing button to second stop.

Tip ejection Remove used tips by pressing ejector button. Change tips as required.

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Measuring unknown volumes (825/835 models only)

The fine volume resolution of the micrometric mechanism allows for precise liquid measurement by aspiration. Remember set volume before starting, dip pipette tip into liquid and turn adjustment button counter-clockwise until unknown quantity is completely aspirated. Measured volume consists in read volume minus initial set volume.

Example: Volume before aspiration: 10.0 µL. Volume read-out after aspiration: 64.7 µL. Liquid aspirated: 64.7-10.0 = 54.7 µL.

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MAINTENANCE

The models were designed to sustain long term operation with minimal maintenance and service requirements. However it is recommended to control instrument performance in accordance with laboratory internal procedures (SOP, GLP, etc) or at least once a year. Defect parts must be replaced, or instrument returned to dealer for repair. Order original spare parts from authorised dealers only.

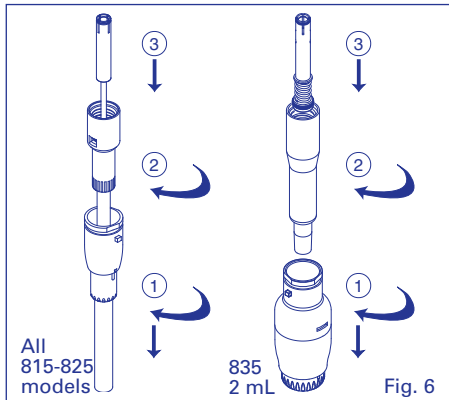
Tightness control

A tight air chamber is paramount for the instrument to provide accurate and precise results. Leaks can be detected by observing a droplet forming and falling from the extremity of the tips or unequal liquid level in the tips of a multichannel model. Special attention must be paid to the following points in case of leakage:

- Matching of tips and nozzle
- Damaged nozzle
- Worn-out PTFE sleeve and/or O-ring

Cleaning

Elements of the volumetric module (lower assembly) can be cleaned with alcohol, or soaked in appropriate decontamination or disinfecting solutions. This operation is particularly important if liquid accidentally entered the pipette. Rinse and dry instrument before reassembling. If ultra-sonic bath is used to remove sticking residues, slightly grease O-ring (see below) and/or plunger before reassembling.

Disassembling volumetric module and O-ring change, 815/825/835 - 2 mL models (Fig. 6)

Removing volumetric module (lower assembly):

1. Press ejector button (2), turn part (20) or (24) for 2 mL to the left and pull.
2. Unscrew barrel (18).
3. Press plunger button in full then unclip (pull) plunger.

O-ring and PTFE sleeve replacement (815/825):

- After removing volumetric module, take barrel (18), press both clips of ring (14) with a pipette tip or a pointed tool.
- Remove ring from barrel, then remove spring (15) to access O-ring/PTFE sleeve assembly.
- Change parts if needed. Clean plunger (13) to eliminate any grease residue.

Note: To prevent damages when reassembling, first place PTFE sleeve (16) on plunger, then O-ring (17) on sleeve.

- Slightly grease plunger, sleeve and O-ring and reassemble volumetric module.

Note: Tightness parts are not available separately on 2, 10, 10Y and 20 µL models. In case of tightness problem, the whole barrel must be changed.

O-ring replacement (835 - 2 mL):

- Unscrew plunger rod (21) and remove washers (22, 23) and spring (15).
- Change parts if needed. Uniformly grease O-ring (17) and barrel (18).

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Disassembling volumetric module and O-ring change, 835 5 and 10 mL models (Fig. 7)

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1. Slightly turn ejector nut (20) to unclip from ejector cap (24).
2. Unscrew barrel (18), press push button in full then pull barrel to unclip.

O-ring replacement:

3. Slightly and simultaneously press both barrel clips (18) with fingers to separate from bonnet (25).
- Pull out plunger assembly.
- Unscrew plunger rod (21) and remove washers (22, 23) and spring (15).
- Change parts if needed. Uniformly grease O-ring (17) and barrel (18).
- Reassemble plunger assembly, barrel and bonnet, screw barrel on pipette, then clip plunger rod by pressing in full plunger button (B).

To use both Pasteur pipettes and PP tips, replace regular barrel by Pasteur pipette adapter nozzle (1.835.633).

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1. Turn ejector nut clockwise till its lowest position.
2. Press ejector button (2), turn ejector nut to the left, unclip ejector ring and release ejector button.
3. Hold volumetric module firmly with one hand and, while pulling down casing, rotate slowly to unscrew.
4. Press plunger button then unclip plunger rod (28).

Barrel change (Fig. 9 to 11):

Note: Tightness O-ring cannot be removed from barrel. Change barrel if tightness is deficient.

1. Press with pointed tool on both clips of cover (27) and remove casing (31).
2. Press bottom clips a) of barrel holder and separate.
3. Pull out barrel assembly without bending the plunger head.

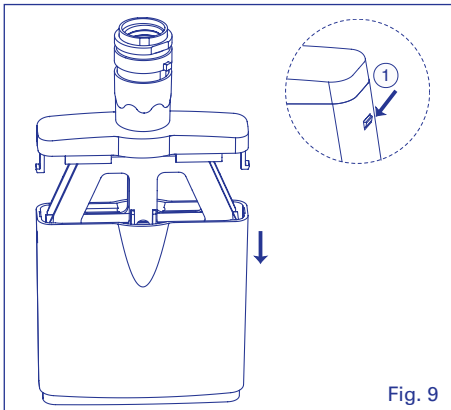
Fig. 9

Reassembling volumetric module, 855 model:

- Apply thin, regular grease layer on all surface length of plunger.
- Separate both plate of barrel holder.
- Introduce barrel (30) on plunger (29) without bending it.
- Clip barrel plate, all barrels must be sitting properly and aligned.
- Introduce casing (31) and clip on cover (27).

Note: Markings on casing (volume) and on cover should not appear simultaneously.

- Hold casing with fingers pulling against barrel extremities.
- Slowly screw volumetric module.
- Press ejector button (2) and introduce tooth of ejector ring in recess.
- Turn counter clockwise to click in ejector rod (11) and release ejector button.
- Press firmly plunger button until plunger rod (28) clips in.
- Control proper functioning of plunger and ejector buttons.
- Set ejector height to desired position.

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

TROUBLE SHOOTING

| Observation | Possible cause | Action |
|--|---|---|
| Tips not fitting the nozzle | Inadequate position of tip ejector Inadequate tips | Adjust position of tip ejector Use original or other compatible tips |
| Pasteur pipette does not hold or too «tight» | Pasteur not compatible | Check/change O-rings Grease slightly O-rings |
| Plunger sticking or dragging | Dirty plunger | Disassemble and clean plunger (835, 855 grease as well) |
| Liquid not aspirated | Dry residues block pipette nozzle Volumetric module not properly assembled Dirty or wet nozzle filter on 835 models | Disassemble and clean Reassemble according to instructions Change nozzle filter |
| Uneven liquid level in tips of multichannel model | Lack of tightness | Check tip fitting Change for damaged nozzle or tightness parts |
| Poor instrument performance | Lack of tightness Inadequate tips Dirty or wet nozzle filter on 835 models Instrument out of calibration Instrument used with viscous or volatile liquids or at particular temperatures | Check for damaged nozzle or tightness parts Check tip compatibility and fitting Pasteur not compatible, check tightness with adapter Change nozzle filter Perform new calibration Perform new calibration with liquid or at temperature used |

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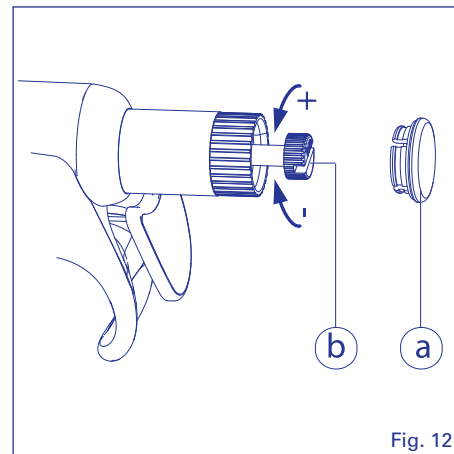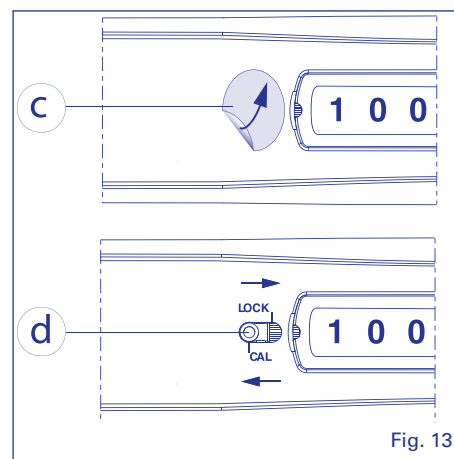Sterilisation

Instrument has been designed for repeated sterilisation in the autoclave at 121°C (20 minutes), fully assembled. Before autoclaving, remove nozzle filter on 835 models. Place instrument horizontally in the autoclave, avoiding any direct contact with metal.

Allow instrument to dry and cool down before use. Insert new nozzle filter on 835 models. Check tightness and accuracy regularly but at least after 50 autoclaving cycles. Check volumetric module, tighten barrel if loose. Repeated autoclaving may affect material colours and free rotation of *smartie* colour caps. Correct autoclaving and resulting sterility are the responsibility of the user.

CALIBRATION

Each Acura® *manual* has been factory calibrated and individually controlled. The QC certificate accompanying the instrument includes control data, serial number and operator identification. Calibration is carried out gravimetrically with aqua-bidest. at constant temperature ($\pm 0.5^\circ\text{C}$) comprised between 20 and 25 °C according to ISO 8655.

In-lab calibration

The *swift-set* calibration system* with integrated key and locking mechanism makes tedious procedures obsolete. If performance results are no longer within recommended values, for instance after QC check or replacing parts, or if changes occur in physical parameters (liquid density, temperature), recalibration is easily performed using an analytical balance. Ten measurements should be first performed on small volume after checking seal tightness and perfect working condition of the instrument.

Adjustable volume models (Fig. 12 and 13)

- Remove *smartie* cap (a) on plunger button and pull out calibration screw (b).
- Remove "Calibration seal" (c) and move calibration cursor (d) on cal position (use pointed item such as tip or clip).
- Turn calibration screw until average weighed volume appears in the display window, without holding push button.
- Place calibration cursor on *lock* position. Press fully to the overtravel position. Cover plunger button with *smartie* cap again.
- Perform few plunger movements and check resulting volume.
- Place a new "Calibration seal" over the calibration cursor.

Fixed volume models (Fig. 12)

Same as above, except

- Turn calibration screw (b) clockwise to decrease and counter-clockwise to increase resulting volume.
- Check resulting volume after each fine tuning, until pipette performance are within acceptable values.

WARRANTY

Your Acura® *manual* pipette is guaranteed against any material or manufacturing defects for the time period specified in its QC certificate. Damages due to non-respect of manufacturer's instructions, safety precautions or autoclaving conditions, as well as material colour alteration are excluded from the warranty. Repair and replacement of parts do not extend warranty time. Should regular maintenance not eliminate a detected defect, return the instrument to the dealer from whom it was purchased after obtaining return authorization.

Decontaminate instrument prior to returning it and with the corresponding form

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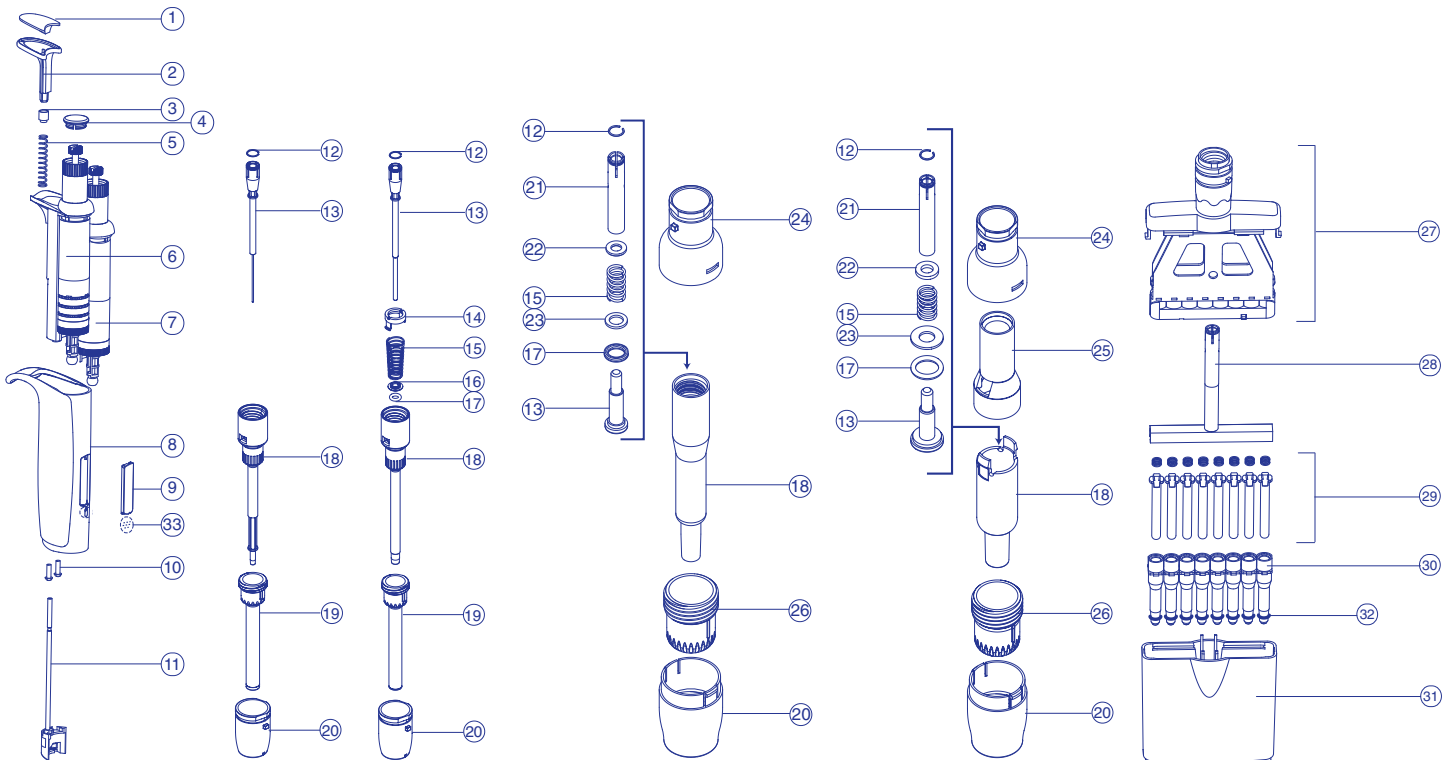
835: 5 and 10 mL

855: all sizes

PERFORMANCE AND ORDERING INFORMATION

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¹⁾ Measured at 0.5 μ L. Indicative data at 0.2 μ L: E< 12%, CV<8%. ²⁾ Measured at 1 μ L. Indicative data at 0.5 μ L: E <+5%, CV<3%.

Acura 835, adjustable volume micropipettes

| Volume mL | Division mL | Inaccuracy (E%) | | | Imprecision (CV%) | | | Tip style | Cat. No. |
|-----------------------|-------------|-----------------|---------------|---------------|-------------------|-----------|-----------|--------------|----------|
| | | Min. vol. | Mid. vol. | Max. vol. | Min. vol. | Mid. vol. | Max. vol. | | |
| 0.2 – 2 | 0.002 | < \pm 1.5 % | < \pm 1.0 % | < \pm 0.5 % | < 0.5 % | < 0.3 % | < 0.2 % | 2 mL | 835.02 |
| 0.2 – 2 ¹⁾ | 0.002 | < \pm 1.5 % | < \pm 1.0 % | < \pm 0.5 % | < 0.5 % | < 0.3 % | < 0.2 % | 2 mL/Pasteur | 835.02PP |
| 0.5 – 5 | 0.01 | < \pm 1.5 % | < \pm 1.1 % | < \pm 0.6 % | < 0.6 % | < 0.5 % | < 0.3 % | 5 mL | 835.05 |
| 0.5 – 5 ¹⁾ | 0.01 | < \pm 1.5 % | < \pm 1.1 % | < \pm 0.6 % | < 0.6 % | < 0.5 % | < 0.3 % | 5 mL/Pasteur | 835.05PP |
| 1 – 10 | 0.01 | < \pm 1.5 % | < \pm 0.7 % | < \pm 0.5 % | < 0.5 % | < 0.3 % | < 0.2 % | 10 mL | 835.10 |

Measurements done with nozzle protection filter. Above values obtained using PP tips.

¹⁾ Model includes Pasteur adapter nozzle.

Acura 855, adjustable volume multichannel pipettes

| Volume μ L | Division μ L | Inaccuracy (E%) | | | Imprecision (CV%) | | | Tip style | Cat. No. |
|-------------------|------------------|-----------------------------|---------------|---------------|-----------------------|-----------|-----------|---------------------|------------|
| | | Min. vol. | Mid. vol. | Max. vol. | Min. vol. | Mid. vol. | Max. vol. | | |
| 8-channel | | | | | | | | | |
| 0.5 – 10 | 0.01 | < \pm 3.5 % ¹⁾ | < \pm 2.5 % | < \pm 1.5 % | < 3.0 % ¹⁾ | < 3.3 % | < 1.0 % | Ultra 10 μ L | 855.08.010 |
| 2 – 20 | 0.02 | < \pm 2.5 % | < \pm 1.8 % | < \pm 1.2 % | < 2.0 % | < 1.4 % | < 0.7 % | Extended 10 μ L | 855.08.020 |
| 5 – 50 | 0.1 | < \pm 1.0 % | < \pm 0.9 % | < \pm 0.8 % | < 1.0 % | < 1.3 % | < 0.4 % | 200 μ L | 855.08.050 |
| 10 – 100 | 0.1 | < \pm 1.0 % | < \pm 0.9 % | < \pm 0.8 % | < 0.9 % | < 0.6 % | < 0.4 % | 200 μ L | 855.08.100 |
| 20-200 | 0.2 | < \pm 0.9 % | < \pm 0.8 % | < \pm 0.7 % | < 0.6 % | < 1.6 % | < 0.3 % | 200 μ L | 855.08.200 |
| 40 – 350 | 0.4 | < \pm 1.0 % | < \pm 0.9 % | < \pm 0.8 % | < 0.6 % | < 1.1 % | < 0.3 % | 350 μ L | 855.08.350 |
| 12-channel | | | | | | | | | |
| 0.5 – 10 | 0.01 | < \pm 3.5 % ¹⁾ | < \pm 2.5 % | < \pm 1.5 % | < 3.0 % ¹⁾ | < 3.3 % | < 1.0 % | Ultra 10 μ L | 855.12.010 |
| 5 – 50 | 0.1 | < \pm 1.0 % | < \pm 0.9 % | < \pm 0.8 % | < 1.0 % | < 1.3 % | < 0.4 % | 200 μ L | 855.12.050 |
| 20-200 | 0.2 | < \pm 0.9 % | < \pm 0.8 % | < \pm 0.7 % | < 0.6 % | < 1.6 % | < 0.3 % | 200 μ L | 855.12.200 |
| 40 – 350 | 0.4 | < \pm 1.0 % | < \pm 0.9 % | < \pm 0.8 % | < 0.6 % | < 1.1 % | < 0.3 % | 350 μ L | 855.12.350 |

¹⁾ Measured at 1 μ L. Performance values obtained with bi-dest. water at constant temperature (\pm 0.5°C) comprised between 20 and 25°C in accordance with ISO 8655.

Acura 835, fixed volume macropipettes

| Volume mL | Inaccuracy (E%) | Imprecision (CV%) | Socorex Cat. No. | Tip style |
|-----------|-----------------|-------------------|------------------|--------------------|
| 2 | <+/- 0.5 % | < 0.2 % | 835.F02 | 2 mL ¹⁾ |
| 2.5 | <+/- 0.8 % | < 0.5 % | 835.F02.5 | 5 mL |
| 5 | <+/- 0.6 % | < 0.3 % | 835.F05 | 5 mL |
| 10 | <+/- 0.5 % | < 0.2 % | 835.F10 | 10 mL |

¹⁾ 5 mL tip up to serial number 17091050

smarties plunger caps

| Colour | Socorex Cat. No. | Colour | Socorex Cat. No. |
|---|------------------|----------|------------------|
| smartie plunger caps, alternative colour selection (6 / pack) | | | |
| White | 825.710 | Rose | 825.719 |
| Grey | 825.715 | Pink | 825.720 |
| Vanilla | 825.716 | Purple | 825.721 |
| Lemon | 825.712 | Ice Blue | 825.722 |
| Yellow | 825.717 | Blue | 825.713 |
| Orange | 825.718 | Mint | 825.723 |
| Red | 825.714 | Green | 825.711 |
| smartie mix assorted colours (14 / pack) | | | 1.825.700 |

ACCESSORIES

ZUBEHÖR

ACCESSOIRES

| Description | Packaging | 7 positions | 6 positions |
|---|-----------|-------------|-------------|
| | | Model 337 | Model 336 |
| Universal work station – fits single channel pipettes of most brands | | | |
| Grey | 1 / pack | 320.337G | 320.336Q |
| Blue | 1 / pack | 320.337B | 320.336B |
| Yellow | 1 / pack | 320.337Y | 320.336Y |
| Green | 1 / pack | 320.337M | 320.336G |
| Pink / Red | 1 / pack | 320.337R | 320.336R |
| Orange | 1 / pack | – | 320.336O |
| White | 1 / pack | – | 320.336W |

| Description | Packaging | Socorex Cat. No. |
|---|------------|------------------|
| Work station 340 for 3 multi- or single-channel pipettes | 1 / pack | 320.340 |
| Shelf pipette holder 332 for 2 instruments (transparent blue) | 1 / pack | 320.332 |
| | 4 / pack | 320.332.4 |
| V-shape, 75 mL, autoclavable PP reservoir | 20 / pack | 330.01 |
| Large capacity with lid, 125 mL, autoclavable PP reservoir | 20 / pack | 330.01L |
| 8-channel, 48 mL, sterile, PS reservoir | 30 / pack | 330.08 |
| | 100 / pack | 330.08XL |
| 12-channel, 48 mL, sterile, PS reservoir | 30 / pack | 330.12 |
| | 100 / pack | 330.12XL |
| 8-channel, 40 mL, sterile, single wrapped PS reservoir | 10 / pack | 330.08.9 |
| 12-channel, 40 mL, sterile, single wrapped PS reservoir | 10 / pack | 330.12.9 |
| Nozzle protection filters, cellulose for 2 and 5mL models | 250 / pack | 322.05 |
| Nozzle protection filters, PP fibres for 10mL model | 100 / pack | 322.10 |
| Pasteur pipette adapter, 2 mL model | 1 / pack | 1.835.631 |
| Pasteur pipette adapter, 5 mL model | 1 / pack | 1.835.633 |
| Protective calibration seal stickers | 25 / bag | 1.835.025 |
| | 100 / bag | 1.835.100 |

QUALITIX® PIPETTE TIPS
QUALITIX® PIPETTENSPITZEN
EMBOUTS QUALITIX®

Superior pipette tips cover a range between 0.2 µl and 10 ml. All models autoclavable at 121 °C. Certified DNase, RNase. Pyrogen and ATP free. Low protein binding tips certified protease free.

| Tips without filter | Description | Vol. (µL) | Racks | Socorex Cat. No. | Bags | Socorex Cat. No. |
|--|--|-----------|--------------|------------------|--------------|------------------|
|  | Natural ultra-microtip | 10 | 12 x 96/rack | 3090.0010R | 1 x 1000/bag | 3090.0010B |
| | Natural ultra-microtip, <i>tipfill</i> rack filling system | 10 | 13 x 96/rack | 3090.0010T | | |
| | Empty rack for <i>tipfill</i> system | 10 | 10/pack | 3070.0210ER | | |
| | Natural ultra-microtip, sterile | 10 | 12 x 96/rack | 3090.0010RS | | |
|  | Natural ultra-microtip, low protein binding | 10 | 12 x 96/rack | 3090.0010PR | | |
| | Natural ultra-microtip, low protein binding, sterile | 10 | 12 x 96/rack | 3090.0010PRS | | |
|  | Natural extended microtip | 10 | 10 x 96/rack | 3020.0010R | 1 x 1000/bag | 3020.0010B |
| | Natural extended microtip, low protein binding | 10 | 10 x 96/rack | 3020.0010PR | | |
| | Natural extended microtip, low protein binding, sterile | 10 | 10 x 96/rack | 3020.0010PRS | | |
|  | Natural gel load tip | 10 | 4 x 204/rack | 302.0010GR | | |
|  | Natural universal tip | 200 | 10 x 96/rack | 3070.0200R | 1 x 1000/bag | 3070.0200B |
| | Natural universal tip, <i>tipfill</i> rack filling system | 200 | 10 x 96/rack | 3070.0200T | | |
| | Empty rack for <i>tipfill</i> system | 200 | 10/pack | 3070.0210ER | | |
| | Natural universal tip, sterile | 200 | 10 x 96/rack | 3070.0200RS | | |
|  | Natural universal tip, low protein binding | 200 | 10 x 96/rack | 3070.0200PR | | |
| | Natural universal tip, low protein binding, sterile | 200 | 10 x 96/rack | 3070.0200PRS | | |
|  | Yellow universal tip | 200 | 10 x 96/rack | 3290.0200R | 1 x 1000/bag | 3290.0200B |
| | Yellow universal tip, sterile | 200 | 10 x 96/rack | 3290.0200RS | | |
|  | Natural gel load tip | 200 | 6 x 96/rack | 3080.0200GR | | |
|  | Natural tip | 350 | 10 x 96/rack | 308.0350R | 1 x 1000/bag | 308.0350B |
|  | Natural universal tip | 1000 | 6 x 96/rack | 3070.1000R | 1 x 1000/bag | 3070.1000B |
| | Natural universal tip, <i>tipfill</i> rack filling system | 1000 | 5 x 96/rack | 3070.1000T | | |
| | Empty rack for <i>tipfill</i> system | 1000 | 10/pack | 3070.1000ER | | |
| | Natural universal tip, sterile | 1000 | 6 x 96/rack | 3070.1000RS | | |
|  | Natural universal tip, low protein binding | 1000 | 6 x 96/rack | 3070.1000PR | | |
| | Natural universal tip, low protein binding, sterile | 1000 | 6 x 96/rack | 3070.1000PRS | | |
|  | Blue universal tip | 1000 | 6 x 96/rack | 3190.1000R | 1 x 1000/bag | 3190.1000B |
| | Blue universal tip, sterile | 1000 | 6 x 96/rack | 3190.1000RS | | |
|  | Natural tip | 1000 | 6 x 96/rack | 3090.1000R | 1 x 1000/bag | 3090.1000B |
|  | Natural macrotip | 2 mL | | | 2 x 250/bag | 312.02 |
|  | Natural macrotip | 5 mL | 2 x 50/rack | 312.05R | 2 x 250/bag | 312.05B |
|  | Natural macrotip | 10 mL | | | 3 x 100/bag | 312.10 |

Efficient protection against aerosols. Include high density PE filter for complete, uncontaminated sample retrieval. Certified DNA, DNase, RNase free. Pyrogen and ATP free

| Filter tips | Description | Vol. (µL) | Racks | Socorex Cat. No. |
|--|--|-----------|--------------|------------------|
|  | Sterile natural ultra-microtip | 10 | 12 x 96/rack | 3090.0010FRS |
|  | Sterile natural ultra-microtip, extended | 10 | 10 x 96/rack | 3020.0010FRS |
|  | Sterile natural tip | 20 | 10 x 96/rack | 3080.0020FRS |
|  | Sterile natural tip | 100 | 10 x 96/rack | 3080.0100FRS |
|  | Sterile natural tip | 200 | 10 x 96/rack | 3070.0200FRS |
|  | Sterile natural tip, extended | 200 | 6 x 96/rack | 3080.0200LFRS |
|  | Sterile natural tip | 300 | 10 x 96/rack | 308.0300FR |
|  | Sterile natural tip | 1000 | 6 x 96/rack | 3070.1000FRS |

B = bag, R = rack, P = low binding, G = gel, L = long, S = sterile, F = filter, T = refill system