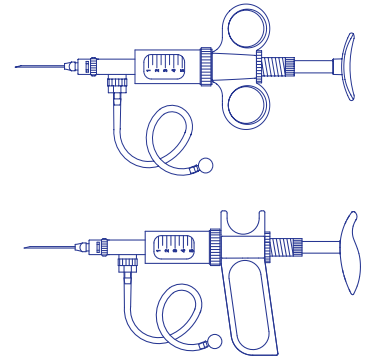
DOSYS™ 162/163/164/172/173/174

LABORATORY SELF-REFILLING SYRINGES LABOR SELBSTFÜLLER-SPRITZEN SERINGUES AUTOMATIQUES DE LABORATOIRE

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SAFETY PRECAUTIONS

- Read the operating instructions carefully before use. Keep available for future reference
- Observe the manufacturer's recommendations
- Follow the regulations on handling of potentially hazardous reagents
- Before each use, check instrument for perfect working condition
- Assure cleanliness when performing cytotoxic sensitive tests
- Prime, rinse and dry the instrument carefully after each use
- Never try to separate blocked parts by using force
- Mind possible hand fatigue during serial dispensing and its medical consequences
- Do not use syringe for purposes other than intended for
- A saline environment may hasten replacement of instruments parts
- Use only original spare parts

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DESCRIPTION

Congratulations on buying a fine quality Swiss product. This self-refilling laboratory syringe possesses a spring-loaded plunger and a 3-way valve and is intended for continuous liquid dosing with a high accuracy and an excellent reproducibility. Good plunger tightness without O-ring, easy volume reading on the interchangeable glass barrel and adjustment mechanism with a locking nut are additional key features of the syringe. All models can be autoclaved fully assembled at 121° C (20 minutes).

The following Dosys™ syringes provide efficient solutions for most repeat dispensing activities including sterile filtration (using accessory kit for enhanced tightness) and dispensing of hot agar or culture media.

Dosys™ basic 162/172 automatic filling and shot by shot dispensing through needle or cannula. Supplied without valve unit. Cover volumes from 0.1 to 10 mL.

Dosys™ classic 163/173 self-refilling and dispensing through integrated 3-way valve. Cover volumes from 0.025 to 20 mL. Most efficient for aqueous, non corrosive reagents. They may be used for small laboratory animal vaccination.

Dosys™ premium 164/174 self-refilling and dispensing through detachable 3-way valve. Cover volumes from 0.1 to 10 mL. Special materials offer enhanced corrosion resistance.

Practical tabletop syringe work station to be ordered separately.

HOLDING THE SYRINGE

The ergonomic shape of the syringe allows a long term repeat dispensing without fatigue.

To operate 162, 163 and 164 models, place forefinger and middlefinger in the two rings of the syringe handle (fig. 1) and action spring-loaded plunger with thumb.

On 172, 173 and 174 models place hand in handle and action plunger with hand palm (fig. 2).

To change position of the handle (41), untighten screw (44), bring handle in the desired position and retighten screw (44).

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OPERATION

The self-refilling syringe is supplied ready for use. If liquid is taken from an opened flask, connect sinker (24) or feed needle (26) to the end of the feed tube (23) and immerse it in the liquid (fig. 3). When liquid is aspirated from a sterile bottle with septum closure, puncture the bottle cap with the vent cannula coupled with a sterile filter. Then connect the (sterilised) feed needle to the end of the feed tube and puncture septum of the bottle.

Volume setting

Tighten locking nut (4) and adjust plunger edge to wanted volume graduation of glass barrel by turning socket (3). Tighten locking nut (4) again in order to prevent any volume modification during activation of the plunger.

Priming

Activate the plunger until the liquid completely fills up the syringe. For efficient priming, set small volume models to their maximal volume. No air bubble must appear in the transparent glass barrel.

Dispensing

Attach a Luer Lock cannula on outlet nozzle and push plunger to dispense liquid. The spring-loaded plunger and three way valve assure an immediate automatic instrument filling. Select internal diameter of the cannula in relation with the distributed volume. A too small internal diameter renders the plunger activation more difficult and reduces the instrument performance.

End of dosing

When dosing is completed, remove the tubing from the feed source and activate the plunger several times in order to eliminate any residual liquid. Disassemble, clean and dry the instrument parts according to the indications in the "maintenance" chapter. **Never store a syringe without prior priming, cleaning and drying.**

Fig. 1

MAINTENANCE

Cleaning

The self-refilling syringe must be carefully cleaned and dried after each use. Therefore, aspirate an appropriate cleaning solution (water or detergent) and rinse thoroughly. Activate the syringe several times with air to remove residual liquid before disassembling.

Disassembling

Remove valve assembly (10) on 164 and 174 models. Hold handle (41) and unscrew casing (8, 9) in order to remove glass barrel (7). Push rod (1) to remove plunger (6). Dry the valve assembly (seat, ball), syringe body, barrel and plunger before reassembling and storing. Using on a regular basis the ABF Premium lubricant after cleaning/drying will protect the plunger and extend the syringe lifespan.

Trouble shooting

A clogged aspiration valve can easily be fixed up by gently detaching the ball from the valve seat. Therefore, introduce a cleaning wire into the intake valve opening and push until the ball moves. If this operation is not possible, immerse the valve in hot water containing an appropriate detergent and operate the plunger. Make sure to rinse the parts thoroughly and dry them completely.

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To prevent a mispositioning of aspiration valve ball, do not release spring loaded plunger rod abruptly while aspirating.

During the distribution, smooth plunger movement and tightness are guaranteed by a thin film of liquid between the plunger and the barrel. Repeat use of the syringe without liquid may damage plunger or barrel. After long activation, some liquid may be found in the back of the plunger and should be removed when dosing is completed.

Only use original spare parts to replace any damaged components of the syringe.

Sterilization

Warning: PVC protection sleeve on glass barrel is not autoclavable. Remove sleeve before autoclaving the instrument.

The laboratory syringe can be autoclaved fully assembled at 121° C, 250° F during 20 minutes (1 ATM). Do not sterilise an instrument that has not previously been primed cleaned and dried. Risk of glass breakage during autoclaving is reduced by slightly untightening the syringe body. Control tightness of all connections before each use.

Correct autoclaving and resulting sterility are the responsibility of the user.

MATERIALS

All parts in contact with the liquid are made of carefully selected, chemically resistant materials:

- Feed tubing: silicone
- Barrel: sodocalcic AR glass
- Plunger: stainless steel, (DIN 316L superior quality on 164 and 174 models)
- Detachable valve body: stainless steel
- Barrel washer: FPM
- Washers for valve; PTFE
- Valve springs and balls: stainless steel
- Sinkers and aspiration cannula: nickel/chrome plated brass

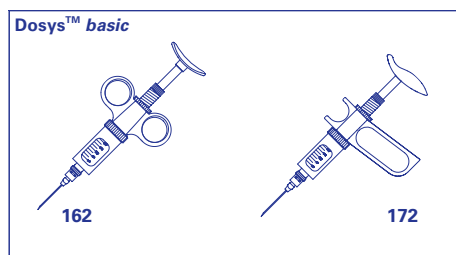
PERFORMANCE AND ORDERING INFORMATION

Package contents

Self-refilling syringe supplied with LL nozzle, silicone tubing, sinker (except 20 mL), aspiration, vent canulas, set of spares for valve (spring, valve balls, washers) cleaning wire and operating instruction. Use vent canula as sinker on 20 mL model.

Superior corrosion resistant Dosys™ premium models 164/174 also include dosing canulas.

Dosys™ basic 162/172 supplied without accessories.

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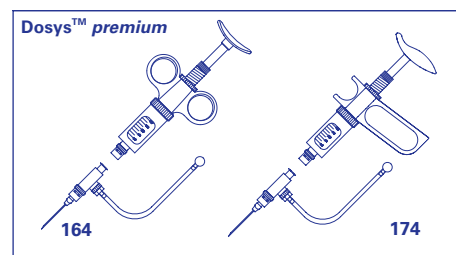
Viscosity Data

Volume mL	Maximal cinematic viscosity* (cst)	Example of liquids
0.3	1800	SAE 50 oil
0.5	2500	Fluid honey
1	1200	Pure Glycerol
2	400 (up to 600) ¹⁾	Glycerol

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Volume mL	Maximal cinematic viscosity* (cst)	Example of liquids
5	150 (up to 200) ¹⁾	SAE 20 oil
10	20	Cream
20	50	Vegetable oil

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Test conditions to determine max. viscosity
Syringe with horizontal feed tubing
Syringe set at maximal volume
Aspiration time < 3 sec

*cinematic viscosity in centistokes (cst). Dynamic viscosity in centipoises (cps) obtained by multiplying the cinematic viscosity by liquid density (kg/dm³)

¹⁾ when using viscous liquid spring (see accessories) ²⁾ at 0.075 mL

ACCESSORIES

Cat. No.	Description	Packaging
Spare glass barrels		
0187003	0.3 mL	6/pk
0187005	0.5 mL	6/pk
018701	1 mL	6/pk
018702	2 mL	6/pk
018705	5 mL	6/pk
018710	10 mL	6/pk
018720	20 mL	3/pk
Spare silicone feed tubings		
187.705	0.3 - 10 mL	1 meter
187.707	20 mL	1 meter
Enhanced plunger tightness set		
1.170.205	For 5 mL model	1 pc
1.170.210	For 10 mL model	1 pc
018705T	Spare barrel for 5 mL set	6/pk
018710T	Spare barrel for 10 mL set	6/pk
Multiple dispensing manifold		
1.170.054	4 channel, LL, SS	1 pc
1.170.058	8 channel, LL, SS	1 pc
Conversion set without valve (not for 164/174)		
1.187.1C	Slant vial holder	1 pc

ZUBEHÖR

Cat. No.	Description	Packaging
Spare parts for syringe valves		
1.187.72003	0.3 mL	1 bag
1.187.72005	0.5 mL	1 bag
1.187.7201	1 mL	1 bag
1.187.7202	2 mL	1 bag
1.187.7205	5 mL	1 bag
1.187.7210	10 mL	1 bag
1.187.7220	20 mL	1 bag
Nylon brush for barrel cleaning		
1.170.301	0.3, 0.5, 1 mL	6/pk
1.170.305	1, 2, 5 mL	6/pk
1.170.320	5, 10, 20 mL	3/pk
1.170.300	set of 2 each	6/pk
Various syringe accessories		
320.170	Work station	1 pc
1.170.950	Detachable valve unit	1 pc
1.170.952	Reverse valve unit	1 pc
187.752	Viscous liquid spring (2 mL)	1 pc
187.753	Viscous liquid spring (5 mL)	1 pc
370.30180	Canula with sprayer nozzle (180 mm)	1 pc
370.30500	Canula with sprayer nozzle (500 mm)	1 pc

ACCESSOIRES

Cat. No.	Description	Packaging
PVC barrel protection sleeves		
1.170.4005	0.3 / 0.5 mL	6/pk
1.170.401	1 mL	6/pk
1.170.402	2 mL	6/pk
1.170.405	5 mL	6/pk
1.170.410	10 mL	6/pk
1.170.420	20 mL	3/pk
PP reservoirs with Rotlauf thread		
1.187.060	60 mL	1 pc
1.187.150	125 mL	1 pc
1.187.250	250 mL	1 pc
Blunt end lab canulas SS, Luer Lock		
370.0840	0.8 x 40 mm	12/pk
370.1250	1.2 x 50 mm	1 pc
370.22100	2.2 x 100 mm	1 pc
370.22150	2.2 x 150 mm	1 pc
Premium lubricant ABF		
1.170.705	70 mL	1 pc
1.170.906	70 mL (pack of 6)	6/pk
1.170.912	70 mL (pack of 12)	12/pk
1.170.755	500 mL	1 pc
1.170.956	500 mL (pack of 6)	6/pk
1.170.962	500 mL (pack of 12)	12/pk

WARRANTY

The self-refilling syringe is guaranteed for a period of one year against any factory or material defect. Cases due to non-respect of safety precautions and manufacturer's instructions are excluded from the warranty. Glass barrel breakage or damages to the plunger due to improper handling or cleaning are not covered by the warranty.

Should regular maintenance not eliminate a detected defect, return instrument to the dealer from whom it was purchased, after having obtained a return authorization.

Decontaminate instrument prior to its return.

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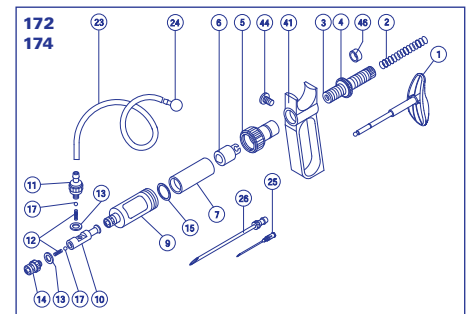
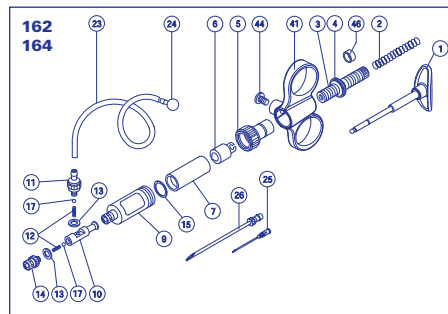
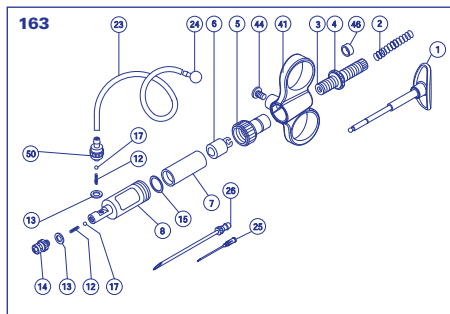
SPARE PARTS

- Piston rod
- Spring
- Socket
- Nut
- Cap
- Piston
- Barrel
- Casing (163/173)
- Casing (164/174, 162/172)
- Valve body (164/174)
- Intake valve (164/174)
- Spring
- Washer
- Nozzle
- Barrel washer
- Valve ball
- Tubing
- Sinker (except 20 mL)
- Vent cannula
- Feed needle
- Handle
- Screw
- PTFE sleeve
- Intake valve (163/173)

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