PROFILLER™ *electro* 447

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SAFETY PRECAUTIONS

- Read operating instructions carefully. Observe manufacturer's recommendations.
- Refer to and follow regulations about handling of potentially hazardous reagents.
- Wear protective clothing including goggles and gloves.
- Do not allow liquid to enter the instrument.
- Do not put aside Profiller™ *electro* 447 if there is liquid in the pipette
- Should the instrument fail to work to specification, stop pipetting and refer to "Trouble Shooting"
- Before cleaning, disconnect power from the charger
- Do not clean the charger.
- All repairs must be carried out by an authorized dealer.
- Only use original accessories and spare parts.
- Do not puncture, crush, bend or modify the battery in any way.

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DESCRIPTION

Profiler™ *electro* 447 is an easy to use and lightweight electronic pipette controller, which facilitates aspiration and distribution of a variety of reagents. The cordless instrument is compatible with glass or plastic pipettes with nominal volumes ranging 0.5 mL to 100 mL or above. The triggers have been designed to eliminate pressure points. Speed control of intake and delivery is therefore made easier.

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APPLICATION LIMITS

- This product should be used only for its intended purpose.
- Profiller™ *electro* 447 is not autoclavable, except silicone pipette holder (G) and nose cone (F).
- Instrument should not be used at temperatures below 10°C and above 35°C.
- Do not use with liquids incompatible with silicone, PP, EPDM and POM.
- Never use with highly flammable liquid, in an atmosphere with danger of explosion, with particular liquid with flash point below 0°C or with acid concentration above 1 mol.
- When instrument is not in use, store on its bench or wall stand in a dry place between -20 °C and +45 °C
- Charger is only to be used indoors and protected against moisture. Only use with manufacturer's original charger.
- Prolonged UV exposure may cause part discoloration. Minimal UV source distance > 50cm.

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Charging

The Profiller™ *electro* 447 operates on Li-ion rechargeable battery.

- Plug the jack directly into the base of the Profiller™ *electro* or into the bench stand (320.948 for Profiller >B00121001).
- Plug the charger after fitting the correct wall adapter (Fig. 1) into a suitable power outlet. Battery charging is indicated with three light bars (Fig. 2).
- Allow unit to charge for 3 to 4 hours from empty to fully charged.
- Fully charged instrument allows for 6h of continuous use.
- Once fully charged, the charging circuit disconnects automatically.
- Batteries service life is approximately 500 charging cycles if used correctly.

Pipetting can also take place while the unit is connected to the charger.

Replacing the battery

If necessary, battery can be replaced easily by user, without any tools (Fig. 1):

1. Take off battery cover and unplug battery cable from socket.
2. Remove battery unit from battery compartment.
3. Slide in new battery unit.
4. Plug in battery cable to socket and close battery cover.

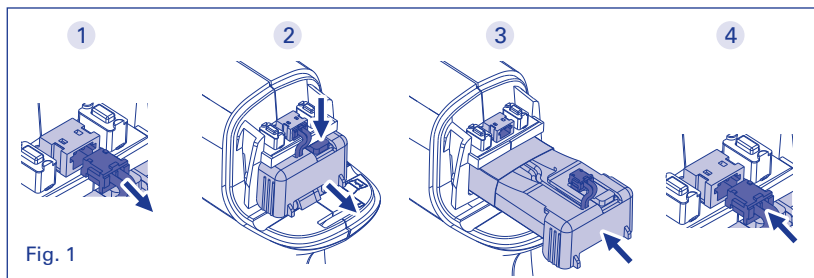
Fig. 1

Notes:

- Before using the instrument for the first time, it is necessary to fully charge the batteries.
- Avoid short charging cycles and disconnect the stand from charger during work when charging is not needed.
- If the indicator shows low battery level during work, connect to the charger to continue working.
- Do not leave the instrument discharged for a long period of time.
- Profiller™ *electro* is not to be disposed with domestic waste. Battery must be disposed in accordance with national regulations.

OPERATION

Speed and mode selection (Fig. 2)

The Profiller™ *electro* 447 has two speeds, HIGH ① and LOW ② set with the SPEED button located on top of the instrument.

It is recommended to set the low speed when working with volumes up to 5 mL.

The MODE button allows two operating modes:

- Button position: ④ Aspirate at selected speed and dispense by gravity (to deliver pipette).
- Button position: ⑤ Aspirate and dispense at selected speed (blow-out pipette).

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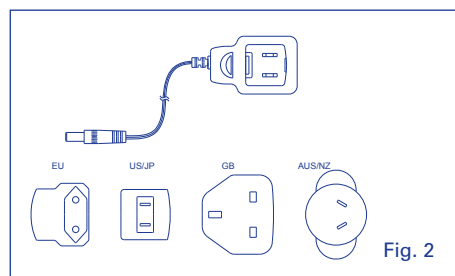
Pipette fitting

Carefully insert glass or plastic pipette into the autoclavable silicone holder of the nose cone without using force.

Aspiration

Hold the instrument in a vertical position with the pipette tip down. Select the speed and operating mode. Place tip of the pipette into the liquid to be aspirated. Press upper trigger (A). The harder the pipetting trigger is pressed the faster the aspiration.

Fill the pipette until the meniscus level has reached the required mark.

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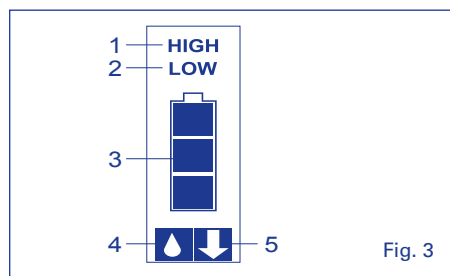
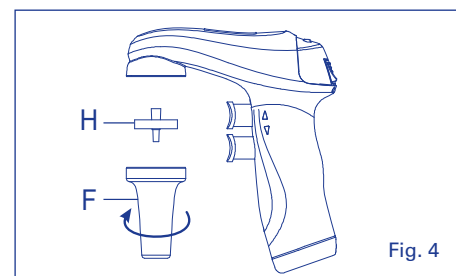
Dispensing

Place pipette against wall of the receiving vessel and press the dispense trigger (B).

In "Blow out" mode: Press the lower trigger (B). The harder the pipetting trigger is pressed the faster the dispensing

In "To deliver" mode: Press the lower trigger (B). Liquid run out by gravity. Wait for the required period of time indicated on the pipette (i.e. EX +10s) for dispensing to be completed.

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Removing the pipette

Hold the pipette and rotate gently while pulling out from the silicone holder.

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TIGHTNESS

Take care of the instrument to keep it in perfect working condition. To test tightness, attach a pipette and fill it at least to half of its nominal capacity. Drops falling from the pipette tip indicate a leak. Unscrew nose cone (F), check silicone pipette holder (G) membrane filter (H) and connector gasket (J). Reassemble correctly.

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MAINTENANCE

Before cleaning disconnect power from charger. Do not clean the charger.

Do not allow liquid to enter the instrument. Hydrophobic protection filter (H) could clog and suction be impaired. To replace filter (Fig 3), remove pipette, unscrew nose cone (F), pull out and replace membrane filter (H). Reassemble unit. Clean instrument with a damp cloth or with most laboratory disinfectants.

Profiller™ *electro* is not autoclavable, except the silicone pipette holder (G) and the nose piece (F) 121°C, 20 minutes.

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Trouble Shooting

The Profiller™ *electro* 447 is virtually service-free. If the recommendations do not resolve the problem, please contact your authorized dealer. Do not attempt to open the instrument.

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TROUBLE	POSSIBLE CAUSE	REMEDY
Aspiration capacity is too slow	Filter (H) blocked, dirty or damaged Batteries (L) low The pipette holder (G) or connector gasket (J) are damaged	Replace hydrophobic filter (0.45 µm or 0.2 µm) Recharge batteries Replace damaged parts
Pipette drips or pipette falls off	Pipette not properly positioned in silicone pipette holder (G) Pipette holder (G) or pipette damaged	Verify pipette mounting Replace damaged parts

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WARRANTY

The Profiller™ *electro* 447 is guaranteed for a one-year period against any factory or material defects. Damages due to non-respect of manufacturer's instructions are excluded from the warranty. Return the instrument to the dealer from whom it was purchased, after obtaining return authorisation, should regular maintenance not eliminate a detected defect.

Decontaminate instrument prior to returning it.

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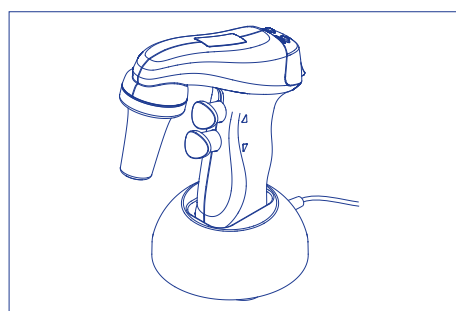
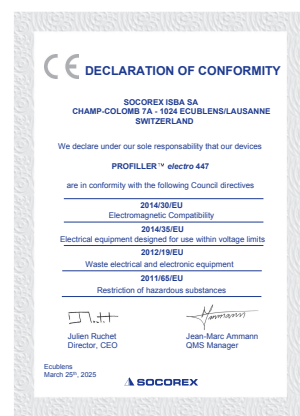
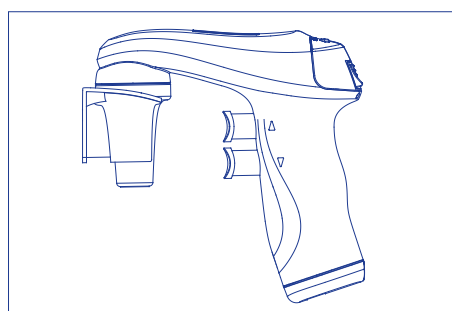
Wall bracket Wandhalterung Support mural

ORDERING INFORMATION

The Profiller™ *electro* 447 is supplied complete with charger, four adaptors (EU, US/JP, GB, AUS/NZ), bench stand and wall bracket, two spare hydrophobic filters (0.45 µm et 0.20 µm) and Quick Guide.

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