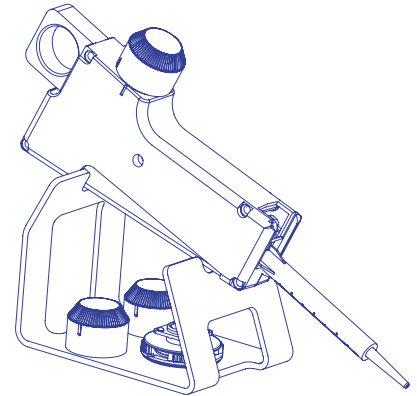
STEPPER™ 416 / ECOSTEP™ SYRINGES

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SAFETY PRECAUTIONS

- Read operating instructions carefully; keep available for future reference
- Observe manufacturer's instructions and guidelines for operation and maintenance
- Refer to and follow regulations about handling of potentially hazardous reagents
- Before use, check instrument for good working conditions
- Use only original Ecostep syringes
- Use water or alcohol to clean instrument (do not use solvents such as acetone)
- Mind possible hand fatigue and its medical consequences
- This product should be used only for its intended purpose

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DESCRIPTION

The Stepper™ 416 is a unique trigger action repeater pipette for fast liquid dispensing without thumb fatigue. Careful design provides for comfort and trouble free dispensing of up to 73 samples without refilling. Sample volume and number of shots are clearly displayed on setting knob. Only three sizes of single use Ecostep™ syringes cover from 10 µL to 5 mL. They are the only part in direct contact with the reagent which greatly reduces risks of carry over or cross contaminations. The Ecostep™ syringes are available bulk or sterile, single wrapped. An automatic locking mechanism prevents dispensing errors when the syringe is almost empty. Performance is better that $\pm 1\%$ ($\pm 2\%$ for volumes under 20 µL).

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Three settings knobs of different colours (one for each syringe size) are supplied with the Stepper. The box includes samples of the Ecostep™ syringes.

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OPERATION

Syringe selection

Non steril Ecostep™ syringe supplied in bulk are perfectly adapted for all routine applications.

When sterility or high purity grade are required, for microbiology or molecular biology applications, choose the single wrapped sterilized Ecostep *bioproof*™. Each production batch is tested to be sterile and free of detectable human DNA, DNase, RNase and pyrogens (endotoxins). Detection levels indicated in batch certificate. See ordering information p.4.

Made of PE/PP, the Ecostep™ syringes show overall good chemical resistance. Use the QR-code below for instant access to the chemical resistance chart. Damages due to incompatible reagents are not covered by the warranty.

Inserting the syringe and setting button (Fig. 1, 2, 3)

- 1 Select button (3) and colour matching syringe (5) according to the volume to be delivered.
 - 2 Insert button (3) on shaft as shown in Fig.1 make sure guide (8) enters slot (9) in button. To remove button, just pull out strongly.
 - 3 Slide rod (2) down until it stops.
- Insert syringe (5) into recess (4) until it stops (Fig. 2).
 - Red syringe needs adapter (7) supplied with the instrument to be fitted first (Fig. 3).

Setting the volume (Fig. 4)

- Unlock setting button (3) by pressing it against casing and rotate until selected volume on scale is even with setting mark (6) on casing.
 - Release button (3) and allow to snap back netally into locking position. Together with the delivered volume, the scale shows the number of shots available when syringe has been filled completely.
- The syringe graduation helps evaluate the number of shots available when the syringe is partially filled.

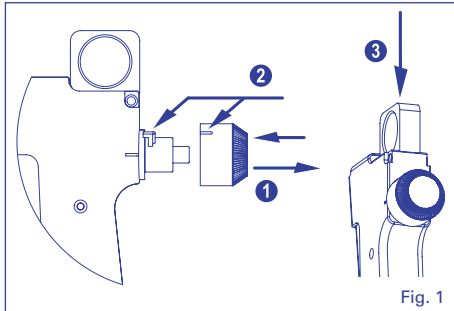
Fig. 1

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DOSING INSTRUCTIONS

Filling the syringe (Fig. 5)

- Immerse syringe nozzle (5) into liquid.
- 1 Fill by **slowly** pulling rod (2) upwards until it stops or, if syringe is to be partially filled, up to the corresponding graduation mark. Small air bubbles on the plunger will not affect precision.
 - 2 Press trigger (1) once until it stops, to position Stepper mechanism on departure line. **Discard first delivery.** When working with expensive reagents, it may be advisable to switch setting button (3) to smallest volume.
- Reposition the setting knob to the desired volume.

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Dispensing (Fig.5)

A pipette tip attached to the syringe will facilitate the access into narrow tubes and increase the delivery precision.

- Place syringe tip against wall of receiving vessel and depress trigger (1) until it stops. Release completely when delivery has been completed
 - To prevent dispensing errors, an automatic device locks the trigger whenever the liquid remaining is less than the set volume. The remaining liquid can be emptied by pushing rod (2) downward.
- Dispensing speed should not exceed 1-2 shots/second.

Removing the Ecostep™ syringe (Fig. 6)

- Push rod (2) downward until it stops.
 - Slide syringe out of recess (4)
- Red syringe adapter can be reused.

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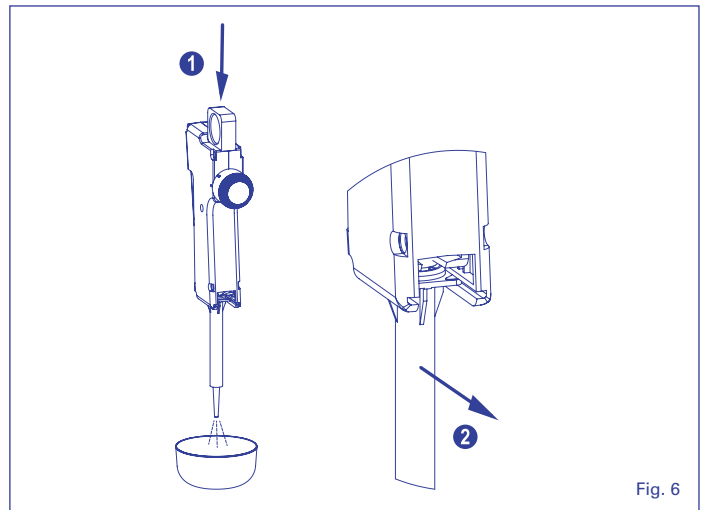
Fig. 6

MAINTENANCE

The Stepper™ 416 does not need any maintenance. Do not open the instrument casing. In case of technical problem, send it to your local dealer for repair after obtaining return authorisation.

Decontaminate instrument prior to returning it.

Calibration

Each Stepper™ is supplied with an individual QC certificate. There is no need and no possibility to calibrate the instrument. Check performance using new, clean Ecostep™ syringes

Sterilization

The Stepper™ and Ecostep™ syringes are not autoclavable. For sterile applications, order sterile, single wrapped Ecostep *bioproof*™ syringes. See page 4 for details.

WARRANTY

The Stepper™ 416 is guaranteed for a period of two years against any factory or material defect. Cases due to non-respect of safety precautions and manufacturers instructions as well as colour alteration are excluded from the warranty. Repair and replacement of parts do not extend the warranty time.

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Chemical compatibility

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**PERFORMANCE DATA
AND ORDERING INFORMATION**

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| Volume | Syringe colour | Syringe capacity | Innaccuracy (E%) | Imprecision (CV%) | Cat. No. |
|---------------|----------------|------------------|--|-------------------|----------|
| 10 – 100 uL | Yellow | 0.75 mL | <+/- 1.5% at 20 uL <+/- 1.0% at 100 uL | < 1.25% < 0.8% | 416.5000 |
| 50 – 500 uL | Blue | 3.75 mL | <+/- 1.0% at 100 uL <+/- 0.5% at 500 uL | < 1.0% < 0.5% | |
| 500 – 5000 uL | Red | 27.75 mL | <+/- 0.8% at 1000 uL <+/- 0.5% at 5000 uL | < 1.2% < 0.4% | |

Performance test in accordance with ISO 8655 method using bi-distilled water at constant temperature (+/- 0.5°C) between 20° and 25°C. Stepper 416 include three colour coded selector buttons, adapter, Ecostep™ syringe samples.

| Ecostep syringe color | Version | Packaging | Cat.No. |
|---|--|---------------|-----------|
| Yellow | Non-sterile | Bulk, 100/box | 316.010 |
| | Sterile <i>bioproof</i> ™ single wrapped | 50 x 1/box | 316.010.9 |
| Blue | Non-sterile | Bulk, 100/box | 316.050 |
| | Sterile <i>bioproof</i> ™ single wrapped | 50 x 1/box | 316.050.9 |
| Red | Non-sterile | Bulk, 100/box | 316.500 |
| | Sterile <i>bioproof</i> ™ single wrapped | 50 x 1/box | 316.500.9 |
| Ecostep selection pack 20 ea. yellow, blue and red syringes | | 60/pk | 316.900 |
| Ecostep <i>bioproof</i> ™ selection pack 20 ea. Sterile, single wrapped yellow, blue and red syringes | | 60 x 1/pk | 316.900.9 |

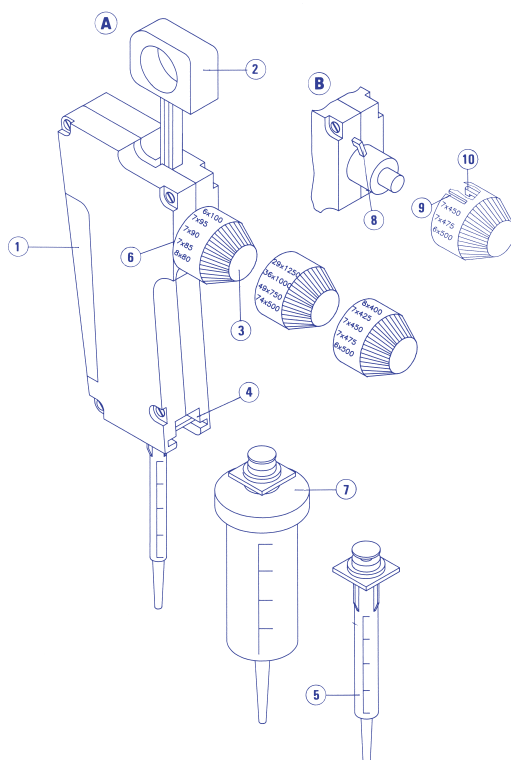
Ecostep *bioproof*™ syringes are certified free of detachable human DNA, DNase, RNase and pyrogen

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